

Instrument Method: MeOH 10%

Stored: 10/8/2018 11:08:36 AM EEST

Method Information

Comments
 Modified User Lauma
 Locked No
 Method Id 2422
 Method Version 2
 Edit User
 Source S/W Info Empower 2 Software Build 2154 DB ID: 741452220

EMD1000 Instrument Setup

Configuration

Mode Operate
 Gas On
 Gas Threshold 100

Function 1

Type	MS Scan	End Time	35.00
Ion Mode	ES+	Scan Time	0.40
Data Format	Centroid	Inter-Scan Delay	0.30
Start Mass	102.00	Cone Voltage	30.00
End Mass	700.00		
Start Time	0.00		

Function 2

Type	MS Scan	End Time	35.00
Ion Mode	ES-	Scan Time	0.40
Data Format	Centroid	Inter-Scan Delay	0.30
Start Mass	102.00	Cone Voltage	30.00
End Mass	700.00		
Start Time	0.00		

W2996 Instrument Setup

Type	W2996	Channel 1 Ratio Wavelength	254.0
Instrument Status	On	Channel 1 Threshold	0.001
Channel Name	W2996	Channel 1 Minimum Ratio	0.001
Start Wavelength	200.0	Channel 1 Maximum Ratio	100.000
End Wavelength	450.0	Channel 1 Filter Type	Hamming
Sampling Rate	1.0	Channel 1 Filter Response	0
Resolution	1.2	Channel 2 Mode	Off
Auto Exposure	Yes	Channel 2 Bandwidth	4.8
Lamp On	Yes	Channel 2 Wavelength	254.0
Interpolate 656 nm	Yes	Channel 2 Offset	0.000
Channel 1 Enable	Off	Channel 2 Ratio Wavelength	254.0
Channel 2 Enable	Off	Channel 2 Threshold	0.001
Exposure Time	15.00	Channel 2 Minimum Ratio	0.001
Channel 1 Mode	Off	Channel 2 Maximum Ratio	100.000
Channel 1 Bandwidth	4.8	Channel 2 Filter Type	Hamming
Channel 1 Wavelength	254.0	Channel 2 Filter Response	0
Channel 1 Offset	0.000	Event Enable	Off

Reported by User: Lauma Laipniece (Lauma)

Report Method: Untitled

Report Method ID 111

Page: 1 of 3

Project Name: Interesterifikacija

Date Printed:

1/28/2019

12:04:03 PM Europe/Riga

Spectral Mode	Data3D	Data Delay Enabled	No
Digital Filter Enabled	No		
Digital Filter Response	1.0		

W2690/5 Instrument Setup

Type	W2690/5
Instrument Status	On
Channel Name	2690/5
Description	
Use Channel Monitor	Off
Monitor Parameter	System Pressure
Stroke Volume	25uL (flow rates <= 0.53 mL/min)
Chart Out	%A
Syringe Draw Rate	Normal
Depth Of Needle	0.0
Degas Mode	On
Pump Mode	Gradient
Flow	0.200
%A	90.0
%B	10.0
%C	0.0
%D	0.0
High Limit	4000.0
Low Limit	0.0
Enable Sample Temp	False
Enable Column Temp	False
Bubble Detect	True
Pre Column Volume	0.0
Sample Temp Target	-1.0
Sample Temp Range	5.0
Spurge A	0.0
Spurge B	0.0
Spurge C	0.0
Spurge D	0.0
Column Temp Target	-1.0
Column Temp Range	5.0
Flow Ramp	2.00
Column choice	No Change
Column Equil Minutes	0.00
Needle Wash	Normal
Switch 1	No Change
Switch 2	No Change
Switch 3	No Change
Switch 4	No Change
Use Events	False
Solvent A	H2O, HCOOH
Solvent B	MeOH
Solvent C	
Solvent D	

W2690/5 Gradient Table

	Time	Flow	%A	%B	%C	%D	Curve
1		0.20	90.0	10.0	0.0	0.0	
2	17.00	0.20	10.0	90.0	0.0	0.0	6
3	28.00	0.20	10.0	90.0	0.0	0.0	6
4	35.00	0.20	90.0	10.0	0.0	0.0	6

Revision History

Version 2 10/8/2018 11:08:36 AM EEST User Lauma

Version 1 10/8/2018 11:06:29 AM EEST User Lauma Method (\\Waters-lmlam3w\Dendrimers 2 : 4059) copied into project. (from Full Audit Trail project)

Version 2 6/15/2017 1:58:17 PM EEST User Lauma

Version 1 6/15/2017 9:17:19 AM EEST User Lauma

Method Version Summaries

	Name	Type	Comments
1	MeOH 10%	Instrument	
2	MeOH 10%	Instrument	Method (\\Waters-lmlam3w\Dendrimers 2 : 4059) copied into project. (from Full Audit Trail project)

Method Version Summaries

	Date	Modified User	Locked	Method Id	Method Version	Source S/W Info
1	10/8/2018 11:08:36 AM EEST	Lauma	No	2422	2	Empow er 2 Softw are Build 2154 DB ID: 741452220
2	10/8/2018 11:06:29 AM EEST	Lauma	No	2421	1	Empow er 2 Softw are Build 2154 DB ID: 741452220

Reported by User: Lauma Laipniece (Lauma)

Report Method: Untitled

Report Method ID 111

Page: 3 of 3

Project Name: Interesterifikacija

Date Printed:

1/28/2019

12:04:03 PM Europe/Riga